

LDS-662-03 2205 8225 08.00 Page 1 of 6 1 August 2019



Fully Automatic Switchover Manifold for Gas Cylinders, Brass, 3000 PSIG (AFAM3000B Series- Laboratory Applications)

SPECIFICATION

Fully Automatic Switchover Manifold

The BeaconMedæs Fully Automatic Switchover Manifold (AFAM3000B Series) accommodates multiple cylinders, equally divided into two banks, and provides an uninterrupted supply of gas for the specific application. The manifold is cleaned, tested and prepared for the indicated gas service and constructed following NFPA, ASME B31.3 and CGA guidelines.

Manifold Description

The AFAM3000B Series manifold operates as a fully automatic switchover control system. It monitors cylinder bank pressure electronically, controlling the switchover and eliminating the need to manually reset levers and valves. The easy to read analog gauges show the delivery and individual bank pressure. A series of lights for each bank indicates whether the bank is "in service", "ready for use", or "bank depleted". At a preset pressure, the system automatically changes from the supply bank to the reserve bank without interruption in gas supply. AFAM3000B Series manifolds operate on 24 VAC power (power transformer provided by BeaconMedæs), but will continue to function during a power failure (without switchover capabilities).

Operation and Design

The AFAM3000B Series consists of a manifold box, an alarm/control module and two supply bank headers (one service and one reserve supply). The manifold box is composed of analog cylinder pressure gauges, regulators, pressure relief valves and normally open solenoid valves. The alarm/control module includes indicator lights (green "In Service", yellow "Ready for Use", and red "Bank Depleted"), a buzzer and a silence button. Each supply bank consists of a header bar with flexible hoses, pressure relief valves and cylinder connections.

After initial power-up and with both banks empty, the red light will be illuminated. The bank that is pressurized first will be considered the "in service" bank and is the bank that supplies the piping system. The green "In Service" light (also referred to as the "Service" bank) will light up, while the cylinder bank on stand-by (or "Secondary Supply") will illuminate the yellow "Ready for Use" light once the bank is pressurized.


On the service bank, the gas flows into the manifold box inlet to the bank pressure switch, continuing onto the primary regulator before heading through a solenoid valve, a check valve and the final line regulator. The delivery pressure is controlled by one line regulator and is field adjustable. The gas exits the line regulator and proceeds past the pressure gauge and into the delivery piping.

The gas on the secondary bank flows into the manifold box inlet, then to the bank pressure switch and the other primary regulator until it reaches a solenoid valve. Since this is the secondary bank, the valve is closed, preventing the secondary bank from flowing. Switchover from the "Service" to "Secondary" side is accomplished when the service bank pressure drops below a predetermined point (the switchover pressure is determined by the pressure switches). The control module then signals the secondary bank solenoid to open while closing the service bank solenoid valve at the same time, allowing it to start flowing without any interruption.

After replacing the empty cylinders, open the cylinder valves. The control module will read the cylinder pressure and automatically place the fresh bank of cylinders into secondary reserve (stand by/ready for use) mode. The yellow light will come on indicating that the new bank is ready for use, and the red "Bank Depleted" light will be extinguished. Replacing the empty cylinders is all that is required to reset the switchover manifold.




Standard Configuration- Brass Construction

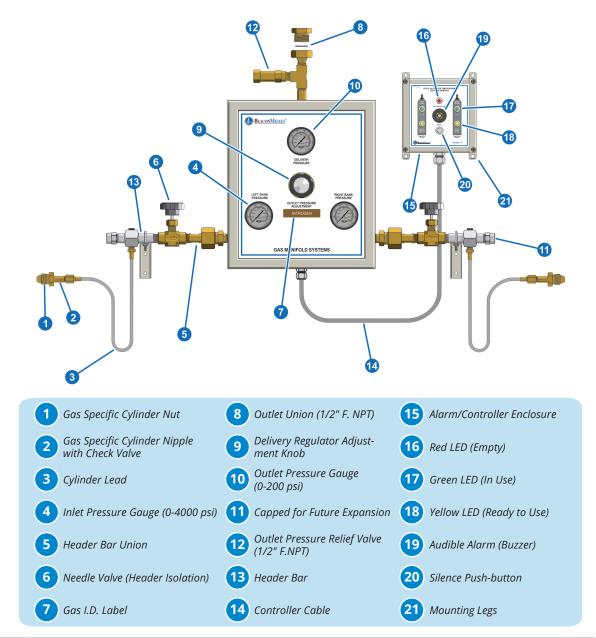


	Enclosure	Steel, Powder Coated, Light Grey	
Materials	Header Bars	Brass Bar Stock, Nickel Plated	
	Tubing	Copper, ASTM B75	
	Fittings	Brass	
	Flexible Hoses	TCH: Teflon Core, Brass Fittings, Stainless Steel Braiding SSH: Stainless Steel Core, Fittings and Braiding	
	Relief Valves	Brass Body, Teflon Seat, Stainless Steel Spring	
	Pressure Reducing Regulators	Brass Body, Stainless Steel Diaphragm, Teflon Seat	
	Pressure Switches	Brass Body, Viton Seals	
	Solenoid Valves	Brass Body, Viton Seals and Seat	
	Optional Vent Valves	Brass Bar Stock, Stainless Steel Diaphragm, PCTFE Seat	



Use & Performance

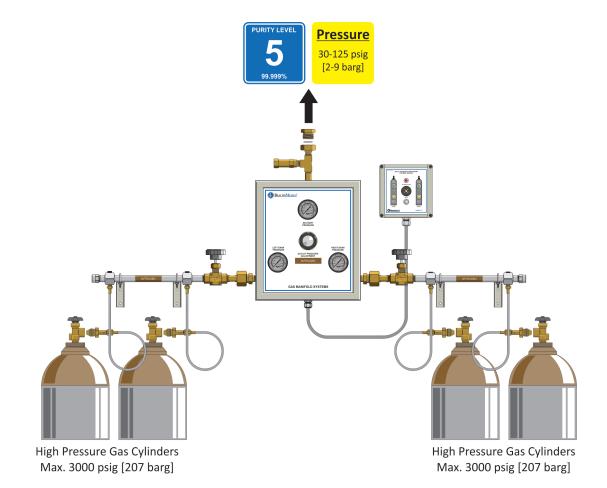

Flow Rate

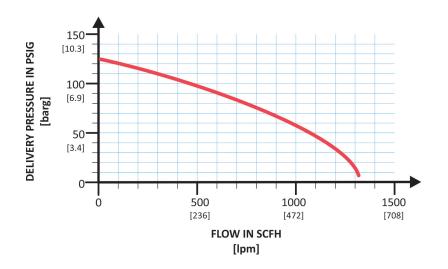



Ordering Information

AFAM3000B - _ - _ - _ - _ - WM - ____ A B C D E F G H

BeaconMedæs AFAM3000 Manifold Model Number Chart						
Variable	Definition	Allowable Value	Description			
Α	Inlet Pressure	3000	3000 PSIG [207 BARG]			
В	Material	В	Brass			
C	Gas	346 580A 580H 350H 580N 590 540	Air- Compressed Argon Helium Hydrogen Nitrogen Air-Industrial Oxygen			
D	No. of Cylinders	2 4 6 8 10	1x1 2x2 3x3 4x4 5x5			
E	Hose	TCH SSH SSHAG	Teflon Core Hose Stainless Steel Hose Stainless Steel Hose with Armour Guard			
F	Configuration	10S 10V	Standard 10" Center Vertical Crossover 10" Center			
G	Mounting	WM	Wall Mount			
Н	Option(s)	VV 3R* IVF	Vent Valves Outside Installation Stainless Steel Tubing for IVF Applications			
(H)	Cylinder Connection	CGA (leave blank) BS DIN NEN	CGA-United States BS341-Great Britain DIN 477-Germany NEN 3268-The Netherlands			

Example: MANIFOLD AFAM, 3000 PSIG INLET, BRASS, HELIUM, 2X2 CYLINDERS, STAINLESS STEEL HOSES, USA *Example Model Number: AFAM3000B-580H-4-SSH-10S-WM*

*3R Option--Outside Installation


AFAM Series Manifolds are designed for indoor installations. If outdoor installation is required, by selecting the "3R" option, all electrical devices (within the manifold and the alarm/controller) will be mounted with NEMA 4X enclosures. Please be aware that even with the 3R option selected, the AFAM enclosure, which is not NEMA 4X, will rust over time due to outdoor exposure as it is made out of steel.



Technical Specifications						
Gas Service	Refer to Part Number Matrix					
Maximum Inlet Pressure	3000 psig [207 barg]					
Delivery Pressure Range	30-125 psig [2-8.6 barg]					
Flow Coefficient	Cv = 0.4, Refer to Flow Chart for details					
Operating Temperature	32°F to 100°F [0°C to 38°]					
Pipeline (Outlet) Relief Valve	200 psig [14 barg]					
Inlet Connections	Gas Specific CGA Fittings					
Manifold Outlet Connection	1/2" F.NPT					
Power	24 VAC (Power Transformer provided by BeaconMedæs)					
Power Transformer	Primary: 120 VAC, Secondary: 24 VAC , 6.25 Amp.					
Maximum Power Consumption	6 Amp					
Electrical Components	All Electrical Components are UL and CSA listed					
Configuration	Normally Open (Supplies Gas When Not Energized)					
Dry Signal Contact	Normally Open (3 Amp. @ 28 VDC/277 VAC when contact is in Close Position					
Cleaning	Cleaned for Oxygen Service as per CGA 4.1					

Remote Alarm Signal Circuitry


The Alarm/Control Box of the AFAM3000B Series Manifold has a dry contact available for remote alarm actuation. It is triggered each time any of the two cylinder banks are empty.

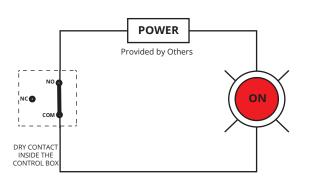

Alarm Condition


When the content inside one of the gas cylinder banks is depleted (low pressure), the dry contact switches from the Normally Closed (NC) position to Normally Open (NO) position. The electrical circuit is closed and the alarm device is actuated.

No Alarm Condition

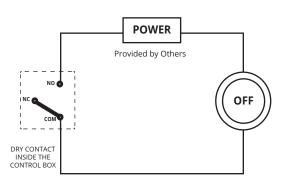

In this situation, both gas cylinder bank pressures are satisfactory (i.e. not empty). The dry contact inside the remote alarm box is in the Normally Closed position. The electrical circuit is open and the alarm device is NOT actuated.



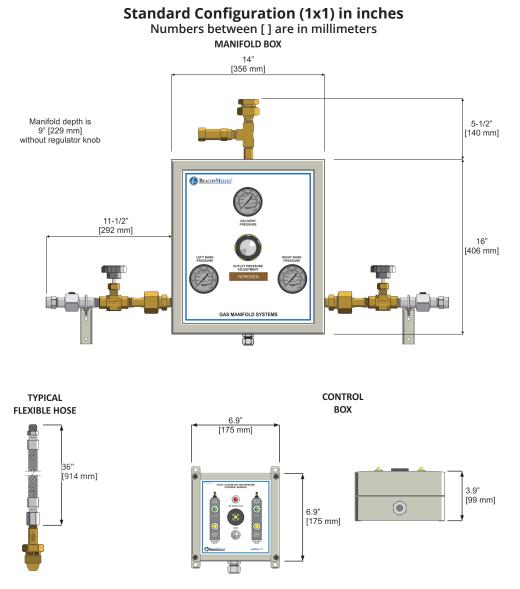

BeaconMedæs AFAM3000 Header Bar Length								
Cylinder	1x1 (10S)	2x2 (10S)	3x3 (10S)	4x4 (10S)				
Configuration								
			5x5 (10V)	7x7 (10V)				
			6x6 (10V)	8x8 (10V)				
Inches	37.50"	57.50"	79.25"	99.50"				
Milimeters	952.5 mm	1460.5 mm	2012.95 mm	2527.3 mm				

BeaconMedæs 1059 Paragon Way, Rock Hill, SC 29730 Phone: +1(803) 817-5600 Fax: +1(803) 817-5750 www.beaconmedaes.com BeaconMedæs Telford Crescent Staveley, Derbyshire, S43 3PF, England Phone: +44(0) 1246-474-242 Fax: +44(0) 1246-472-982 www.beaconmedaes.com